

The Steraffirm Steam Class 6 Emulating Indicators are cycle verification indicators that confirm the presence or absence of specific time and temperature parameters during a cycle, and integrate all the critical parameters of the steam sterilization cycle (temperature, saturated steam and exposure time).



Steraffirm® Steam Class 6 Emulating Indicators

Features

Abrupt and distinctive total color change from yellow to blue/purple

Endpoint color standard is present on each indicator strip

Indicator can be kept as a permanent record

Lot number on each indicator strip and on each box

Conforms to ISO 11140-1 (Class 6)

Benefits

Delivers easy-to-interpret results

Indicates when time and temperature parameters are achieved




Provides documentation of the efficacy of the sterilization cycle

Enhances traceability

Meets critical standards as an indicator for monitoring specific cycle parameters



Emulating Indicators can be placed inside each pack or pouch during sterilization of supplies or equipment in a validated prevacuum cycle. The indicators must be checked by the end user (before using the processed item) to confirm that the specific sterilizing parameters have been attained within the pack. Class 6 emulating indicators can also be placed in test packs and processed in a cycle. The test packs should be opened immediately after the cycle to assess the success of the process. These emulating indicators can be used for routine monitoring of research or pharmaceutical steam sterilization cycles. The Steraffirm Class 6 Emulating Indicators are for life sciences applications only. They are not approved for use in healthcare facilities in the United States.

COLORS ARE AN APPROXIMATION






Steraffirm™ **STEAM** Emulating Indicator

RE ORDER No. PCC012 Class 6
For **STEAM** sterilizers working at 121°C / 250°F for 15 minutes
ACCEPT ONLY IF YELLOW CIRCLE TURNS BLUE/PURPLE



STERIS®


UNUSED









Steraffirm™ **STEAM** Emulating Indicator

RE ORDER No. PCC012 Class 6
For **STEAM** sterilizers working at 121°C / 250°F for 15 minutes
ACCEPT ONLY IF YELLOW CIRCLE TURNS BLUE/PURPLE

STERIS®


TYPICAL PASS





TYPICAL FAIL

Steraffirm® Steam Class 6 Emulating Indicators

Ordering Information

Product	Description	Quantity
PCC012	Steraffirm Steam Class 6 Emulating Indicator 121 °C / 250 °F for 15 minutes	100 / box