

---

## GENERAL DESCRIPTION

CIP Additive™ Detergent Booster and Defoamer is a neutral pH detergent additive for use in boosting the cleaning ability of either alkaline or acidic cleaners. CIP Additive Detergent Booster and Defoamer can be used alone or in combination with other cleaners to supplement surfactancy and to enhance cleaning at temperatures ranging from ambient to 85°C (185°F). CIP Additive Detergent Booster and Defoamer, when used at 43°C (110°F) or higher, can be used in high-energy applications such as recirculating spray or eductor return systems.

---

## FEATURES

Neutral pH detergent additive  
Controlled foaming at higher temperature

---

## BENEFITS

Boosts the cleaning ability of either alkaline or acidic products  
May be used in recirculating systems at temperatures >43°C (110°F) to boost cleaning

---

## PHYSICAL PROPERTIES

Form .....	Colorless to light yellow, clear liquid
Odor.....	None
Specific Gravity (25°C [77°F]).....	1.02, typical
pH (Undiluted) .....	8.57, typical
Solubility .....	Complete
Foam.....	Low at temperatures above 43°C (110°F)
Phosphates .....	None

---

## DIRECTIONS FOR USE

As a detergent additive, the typical use-concentration of the CIP Additive Detergent Booster and Defoamer is 1-4 fl oz/gal (0.8-3.1% v/v). CIP Additive Detergent Booster and Defoamer cannot be titrated. Add in proportion to alkaline or acidic cleaner. Contact STERIS for testing methods for residuals of CIP Additive Detergent Booster and Defoamer.

---

## STORAGE

This product should be stored in an area where it is not exposed to extreme temperatures. Mixing may be needed if stored above 33°C (91°F). Product may freeze. Swirls and precipitate may occur in frozen material, but product will readily go into solution when thawed and mixed.

---

## DISPOSAL

Flush with plenty of water to the sanitary sewer. Dispose of in accordance with local, state, and Federal regulations.

---

## SERVICE

### Sales

Service is one of the most important ways to verify consistent quality of the facility's performance and operation. A tailored service program by STERIS provides effective, trouble-free operations.

## Technical

STERIS is pleased to provide a completely staffed and equipped technical service laboratory capable of performing needed tests and providing both telephone and on-site assistance when needed. More details on how this service can benefit a facility's particular situation can be provided upon request.

## PRECAUTIONS

---

Information concerning human and environmental exposure may be reviewed on the Safety Data Sheet (SDS) for the product. For additional information regarding incidents involving human and environmental exposure, call 1-800-424-9300.

---

### For further information, please contact:



STERIS Corporation  
5960 Heisley Road  
Mentor, OH 44060-1834 • USA  
440-354-2600 • 800-444-9009  
[www.STERISLifeSciences.com](http://www.STERISLifeSciences.com)