



# BIOHAZARD CLEAR BAGS FOR STEAM DECONTAMINATION CYCLES TECHNICAL DATA

---

## GENERAL DESCRIPTION

Biohazard Clear Bags are made of strong polypropylene. They are autoclavable and are labeled with bold visuals signifying decontamination use.

---

## APPLICATION

Disposable, nonsharp items contaminated with biohazardous materials may be placed in Biohazard Clear Bags. Prior to disposal, the filled Biohazard Clear Bags are decontaminated by autoclaving at a temperature and exposure time per the Centers for Disease Control (CDC) guidelines, National Institutes of Health (NIH) guidelines or other appropriate guidelines.

---

## FEATURES

Two sizes available  
Strong material  
Clear plastic

---

## BENEFITS

Accommodate a variety of materials  
Resists tears  
Contents of the bag may be easily verified

---

## TECHNICAL PROPERTIES

**Material:** polypropylene

**Thickness of NY701:** 1.5 mils

**Thickness of NY702:** 2 mils

---

## DIRECTIONS FOR USE

Select the proper size Biohazard Clear Bag. Place the disposable, nonsharp biohazardous items in the bag. When the bag is 3/4 full, seal the bag. Decontaminate with an appropriate steam sterilization cycle.

*NOTE: Standard sterilization cycles are not always adequate for processing highly contaminated materials. Cycle times, temperatures and other conditions should be established for the specific application to ensure sterilization.*

---

## STORAGE CONDITIONS

Biohazard Clear Bags should be kept dry and stored at room temperature.

---

## SERVICE

### Sales

Service is one of the most important ways to verify consistent quality of the facility's performance and operation. A tailored service program by STERIS provides effective, trouble-free operations.

### Technical

STERIS is pleased to provide a completely staffed and equipped technical service laboratory capable of performing needed tests and providing both telephone and on-site assistance when needed. More details on how this service can benefit a facility's particular situation can be provided upon request.

---

## ORDERING INFORMATION

Biohazard Clear Bag Size	Quantity Per Case	Reorder Number
12 x 24" (310 x 610 mm)	200	NY701
24 x 30" (610 x 760 mm)	200	NY702

---

**For further information, please contact:**

**STERIS**<sup>®</sup> STERIS Corporation  
5960 Heisley Road  
Mentor, OH 44060-1834 ■ USA  
440-354-2600 ■ 800-548-4873  
www.steris.com

